

Proteinase K, lyophilized

LOT: See product label

EXPIRY DATE: See product label

ORDERING INFORMATION

CAT.NO.	SIZE	PACKAGE CONTENT
BR1100901	30 mg	30 mg Proteinase K lyophilized
BR1100902	150 mg	5 × 30 mg Proteinase K lyophilized

COMPONENT	COMPOSITION
Proteinase K lyophilized	Lyophilized Proteinase K powder. Add 1.5 ml water to 30 mg Proteinase K.
STORAGE	Store lyophilized Proteinase K at 4°C (until expiry date – see product label) Store aliquots of dissolved Proteinase K at –20°C

FEATURES

- Exceptional purity suitable for the most sensitive applications
- PCR-grade for highest performance
- Stable lyophilizate

APPLICATIONS

- Nonspecific protein degradation
- Nucleic acid purification
- Deactivation of nucleases in enzymatic reactions

Proteinase K, lyophilized

DESCRIPTION

biotechrabbit™ Proteinase K is an active endopeptidase that is effective with native proteins, allowing endogenous RNases and DNases to be inactivated rapidly. The robust enzyme is stable over a wide pH range (4–12.5) and remains fully active for several hours when incubated at pH 6.5–9.5. The exceptional purity of the enzyme ensures that it is ideally suited for preparing PCR templates, as it is free of RNases, DNases and DNA. biotechrabbit Proteinase K is available as an aqueous solution or lyophilized powder.

USEFUL HINTS

Inhibitors

The enzyme is not inactivated by metal ions, chelating agents (such as EDTA), sulfhydryl reagents, or trypsin and chymotrypsin inhibitors.

Activators

The Proteinase K activity is stimulated by denaturing agents (SDS and urea). The presence of SDS can produce a seven-fold increase in Proteinase K activity.

pH Optimum

Proteinase K is stable over a broad pH range (4–12.5). The enzyme retains its full activity for several hours when incubated at a pH between 6.5 and 9.5.

Proteinase K, lyophilized

CERTIFICATE OF ANALYSIS

Quality Control

Protein Purity

The Protein purity is analyzed by SDS polyacrylamide gel electrophoresis.

Absence of nuclease activity

The enzyme is tested for absence of exonuclease, endonuclease and nicking activities.

E. coli DNA contamination assay

The enzyme is tested for absence of *E. coli* genomic DNA.

Quality confirmed by: Head of Quality Control

SAFETY INSTRUCTIONS

For safety instructions please see Safety Data Sheets (SDS)/Sicherheitshinweise finden Sie in den SDS unter:
<http://www.biotechrabbt.com/support/documentation.html>.

USEFUL HINTS

- Visit Applications at www.biotechrabbt.com for more products and product selection guides.
- Most biotechrabbt products are available in custom formulations and bulk amounts.

CONTACT BIOTECHRABBIT

biotechrabbt GmbH
Volmerstr. 9a
12489 Berlin, Germany

info@biotechrabbt.com
support@biotechrabbt.com
www.biotechrabbt.com

Phone: +49 30 555 7821-10
Fax: +49 30 555 7821-99



Legal Disclaimer and Product Use Limitation

Purchase of product does not include a license to perform any patented applications; therefore it is the sole responsibility of users to determine whether they may be required to engage a license agreement depending upon the particular application in which the product is used. This product was developed, manufactured, and sold for in vitro use only. It is not suitable for administration to humans or animals.

Trademarks: biotechrabbt™ (biotechrabbt GmbH).

valid from 24.08.2016