

DNA Blunting Kit

LOT: See product label

EXPIRY DATE: See product label

ORDERING INFORMATION

CAT. NO.	SIZE	PACKAGE CONTENT
BR1101101	100 rxn of 25 µl	125 µl 20× Blunting Enzyme Mix 250 µl 10× Blunting Buffer 25 µl 10 mM dNTP Mix

COMPONENT

COMPOSITION

Blunting Enzyme Mix	20× Blunting Enzyme Mix, in storage buffer containing 50% (v/v) glycerol.
---------------------	---

Blunting Buffer	Optimized 10× Blunting Buffer for rapid blunting reaction.
-----------------	--

10 mM dNTP Mix	dNTP Mix containing 10 mM dATP, dCTP, dGTP, dTTP in water.
----------------	--

STORAGE

-20°C (until expiry date – see product label)

FEATURES

- Rapid blunting reaction
- Excellent performance in blunt-ended ligation
- Ready-to-use blunted product in subsequent ligation

APPLICATIONS

- Blunting and phosphorylation of double-stranded DNA
- Preparation of restriction enzyme digested DNA or sheared DNA for blunt-ended ligation

DNA Blunting Kit

DESCRIPTION

biotechrabbit™ DNA Blunting Kit shows excellent performance in blunting and phosphorylation of double-stranded DNA. The kit is compatible with a wide variety of dsDNA templates such as restriction enzyme digested DNA fragments, plasmid DNA, PCR products with dA-overhang, sheared or nebulized DNA. The kit is optimized for rapid blunting reaction.

PROTOCOL

Blunting reaction

- Scale the reaction volume up as required.
- Prepare the blunting reaction as shown below:

COMPONENT	VOLUME	FINAL CONCENTRATION
10× Blunting Buffer	2.5 µl	1×
Purified DNA	Variable	up to 5 µg
10 mM dNTP Mix	0.25 µl	0.1 mM
20× Blunting Enzyme Mix	1.25 µl	1×
Nuclease free water	Variable	
Total volume	25 µl	

- Mix and centrifuge briefly to collect the liquid in the bottom of the tube.
- Incubate at 25°C (room temperature) for 15 minutes.
- For better efficiency or sheared/nebulized DNA extend the incubation time to 30 min.
- Heat inactivate the enzyme at 70°C for 10 minutes.
- Blunted DNA can be directly used in ligation step (overnight ligation suggested).

DNA Blunting Kit

CERTIFICATE OF ANALYSIS

Quality Control

Functional assay

Purified double digested sticky-end DNA is blunted, ligated and successful ligation is confirmed.

Exonuclease Activity

Linearized lambda/HindIII DNA fragments are incubated with the enzyme in a 50 µl reaction mixture for 4 h at 37°C. No DNA degradation observed.

Endonuclease/Nick Activity

Supercoiled plasmid DNA is incubated with the enzyme in a 50 µl reaction mixture for 4 h at 37°C. No conversion of covalently closed circular DNA to nicked DNA detected.

Contamination with *E. coli* DNA

Absence of *E. coli* genomic DNA is confirmed by qPCR using a sample of the enzyme and specific primers targeting the *E. coli* 16S rRNA gene. No contamination detected.

Quality confirmed by: Head of Quality Control

SAFETY INSTRUCTIONS

For safety instructions please see Safety Data Sheets (SDS)/Sicherheitshinweise finden Sie in den SDS: <http://www.biotechrabbit.com/support/documentation.html>.

USEFUL HINTS

- Visit Applications at www.biotechrabbit.com for more products and product selection guides.
- Most biotechrabbit products are available in custom formulations and bulk amounts.

CONTACT BIOTECHRABBIT

biotechrabbit GmbH

Volmerstr. 9
12489 Berlin,
Germany

info@biotechrabbit.com
support@biotechrabbit.com
www.biotechrabbit.com

Phone: +49 30 555 7821-10
Fax: +49 30 555 7821-99



Legal Disclaimer and Product Use Limitation

Purchase of product does not include a license to perform any patented applications; therefore it is the sole responsibility of users to determine whether they may be required to engage a license agreement depending upon the particular application in which the product is used. This product was developed, manufactured, and sold for in vitro use only. It is not suitable for administration to humans or animals.

Trademarks: *biotechrabbit™* (biotechrabbit GmbH).

valid from 09.04.2019