

RNase Inhibitor, 40 U/ μ l

LOT: See product label

EXPIRY DATE: See product label

ORDERING INFORMATION

CAT. NO.	SIZE	PACKAGE CONTENT
BR0400901	2500 U (125 rxn)	62.5 μ l RNase Inhibitor
BR0400902	10000 U (500 rxn)	4 \times 62.5 μ l RNase Inhibitor

COMPONENT

COMPOSITION

RNase Inhibitor

RNase Inhibitor, 40 U/ μ l, in storage buffer containing 50% (v/v) glycerol.

STORAGE

-20°C (until expiry date – see product label)

FEATURES

- Exceptionally pure proprietary Ribonuclease Inhibitor for demanding RNA applications
- Active under variety of reaction conditions used for work with RNA
- Prevention of RNA from degradation by a wide range of RNases

APPLICATIONS

- In vitro transcription/translation
- cDNA synthesis
- RNA purification and storage

RNase Inhibitor, 40 U/ μ l

DESCRIPTION

biotechrabbit™ RNase Inhibitor is an acidic protein that is a potent inhibitor of a wide spectrum of ribonucleases. The RNase Inhibitor helps to prevent RNA degradation in applications like cDNA synthesis, RT-PCR, in vitro transcription/ translation reactions or RNA purification. The enzyme is purified from a recombinant *E. coli* strain carrying the RNase Inhibitor gene.

PROTOCOL

General guidelines

- RNase Inhibitor can be used in all common reactions performed with RNA, it is active in all common buffers.
- The recommended final concentration for RNA protection is about 1-2 units of RNase Inhibitor for every 1 μ l of the reaction mixture.
- Optimal temperature for Ribonuclease Inhibitor activity is 37°C
- Inactivation conditions: 65°C for 20 minutes

Prevention of cDNA synthesis reaction contamination

RNase contamination is an exceptional concern when working with RNA. RNase A, providing most threat to RNA integrity, is a highly stable contaminant of any laboratory. To prevent RNA from degradation and to minimize possibility of contamination during cDNA synthesis; follow the guidelines below:

- Use separate clean areas for preparation of the samples and the reaction mixture.
- DEPC-treat all tubes and pipette tips or use certified nuclease-free labware with aerosol filters.
- Wear fresh gloves when handling RNA and all reagents.
- Always assess the integrity of RNA prior to cDNA synthesis in denaturing agarose gel electrophoresis.
- Use RNase free water and other reagents.
- To prevent RNA from degradation, add Ribonuclease inhibitor (optional) in to the cDNA synthesis reaction (20 units for 20 μ l reaction).

Typical cDNA synthesis reaction set up

- Thaw on ice and mix very well all reagents.
- Assemble and keep all reactions on ice.
- To use time and reagents effectively, always prepare master mix for multiple reactions. For a master mix volume, always calculate the number reactions that you need plus one additional.
- Combine the following in an RNase-free reaction tube:

COMPONENT	VOLUME	FINAL CONCENTRATION
dNTP Mix (10 mM each dNTP)	2 μl	1 mM (each dNTP)
RNase Inhibitor, 40 U/μl	0.5 μl	1 U/μl
<i>Oligo (dT)₁₂₋₁₈ (10 μM) – or</i>	<i>0.5 μl</i>	<i>0.25 μM</i>
<i>Hexamer Primer (25 μM) – or</i>	<i>1 μl</i>	<i>1.25 μM</i>
<i>Gene Specific Primer (10 μM)</i>	<i>0.5 μl</i>	<i>0.25 μM</i>
RNA Template	0.1–1 μg total RNA or 50–500 ng mRNA (polyA)	
Reverse Transcriptase, 200 U/μl (i.e. BR0400301)	1 μl	10 U/μl
5× RT Buffer	4 μl	1×
RNase-free water (i.e. BR1900301)	Variable	
Total volume	20 μl	

- Mix and collect the drops by centrifuging briefly.
- When using
 - Hexamer Primer, incubate 10 minutes at 30°C followed by 50–55°C for 20–60 minutes
 - Oligo (dT) or gene-specific Primer incubate at 50–55°C for 20–60 minutes.
- Inactivate enzymes at 99°C for 5 minutes.
- Collect the drops by spinning briefly.
- Store products at –20°C or proceed to next step, like PCR or qPCR.
- *Use maximum 10 μl of the cDNA synthesis reaction mix for PCR in 50 μl volume.*

RNase Inhibitor, 40 U/ μ l

CERTIFICATE OF ANALYSIS

Unit Definition

One unit is defined as the amount of enzyme required to inhibit 50% of the activity of 5 ng RNase A (hydrolysis of cyclic cytidine-monophosphoric).

Quality Control

Protein Purity

The Protein purity is analyzed by SDS polyacrylamide gel electrophoresis.

RNase Assay

A sample of the enzyme was incubated with a RNA template. RNase activity was not observed after agarose gel electrophoresis.

Quality confirmed by: Head of Quality Control

SAFETY INSTRUCTIONS

For safety instructions please see Safety Data Sheets (SDS)/Sicherheitshinweise finden Sie in den SDS unter: <http://www.biotechrabbit.com/support/documentation.html>.

USEFUL HINTS

- Visit Applications at www.biotechrabbit.com for more products and product selection guides.
- Most biotechrabbit products are available in custom formulations and bulk amounts.

CONTACT BIOTECHRABBIT

biotechrabbit GmbH
Volmerstr. 9a
12489 Berlin, Germany

info@biotechrabbit.com
support@biotechrabbit.com
www.biotechrabbit.com

Phone: +49 30 555 7821-10
Fax: +49 30 555 7821-99



Legal Disclaimer and Product Use Limitation

Purchase of product does not include a license to perform any patented applications; therefore it is the sole responsibility of users to determine whether they may be required to engage a license agreement depending upon the particular application in which the product is used. This product was developed, manufactured, and sold for in vitro use only. It is not suitable for administration to humans or animals.

Trademarks: biotechrabbit™ (biotechrabbit GmbH).

valid from 24.08.2016